

Pyros[®] eXpress

Endotoxin & Glucan Analysis Software

Release Notes v1.3.74

Pyros[®] eXpress Software supports all of the quantitative endotoxin and glucan detection assays from Associates of Cape Cod, Inc. (ACC), and allows users to quickly and efficiently test in a Quality Control environment. The Pyros[®] eXpress Software provides greater flexibility and versatility in the laboratory allowing you to work smarter and faster while maintaining regulatory compliance. Pyros[®] eXpress Software meets 21CFR Part 11 technical requirements for electronic records, signatures, audit trails as well as US and EU data integrity expectations.

Version 1.3.74

Fixed Issues

- Resolved a security vulnerability that could allow authenticated users with local access to elevate privileges within the application.

CVSS Vector: CVSS:4.0/AV:L/AC:L/AT:N/PR:L/UI:N/VC:H/VI:H/VA:N/SC:N/SI:N/SA:N
CVSS v4.0 Score: 8.4

- Strengthened authentication controls by increasing the minimum password length requirement for application-managed accounts.
- Improved error handling and communication robustness when interfacing with the Pyros Kinetix[®] Flex reader.
- Implemented additional updates to enhance overall system stability, reliability, and security.
- Removed MVC from the calculator when using units of EU/device